


LT-D1201C

1,1':4'1"-terphenyl

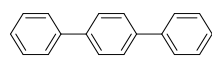


Formula	: C <sub>18</sub> H <sub>14</sub>	Dye Laser Characteristics				
MW	: 230.31 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 92-94-4	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 276 nm(cyclohexane)	Pump source (nm)				
Fluorescence (λ-max)	: 337 nm(cyclohexane)	XeCl(308)	340	334-345	0.46	cyclohexane
		KrF(248)	338	326-359	0.23	cyclohexane

Reference : 1.Appl.Phys.,19, 35(1979)  
2.Optics Commun., 26(3), 396(1978)

LT-D1202C

1,1':4',1":4",1""-quaterphenyl

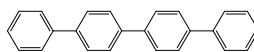


Formula	: C <sub>24</sub> H <sub>18</sub>	Dye Laser Characteristics				
MW	: 306.41 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 135-70-6	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 294 nm(cyclohexane)	Pump source (nm)				
Fluorescence (λ-max)	: 365 nm(cyclohexane)	N2(337)	374	362-390	2.1	toluene

Reference : 1.Nature, 225, 544(1970)

LT-D1203C


3,5,3""',5""'-tetra-*t*-butyl-*p*-quinquephenyl

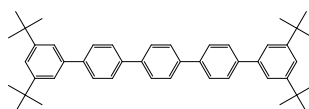

Formula	: C <sub>46</sub> H <sub>54</sub>	Dye Laser Characteristics				
MW	: 606.93 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 89703-16-2	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 311 nm(p-dioxane)	Pump source (nm)				
Fluorescence (λ-max)	: 390 nm(p-dioxane)	N2(337)	387	372-411	0.52	

LT-D1204C

2-methyl-5-*t*-butyl-*p*-quaterphenyl

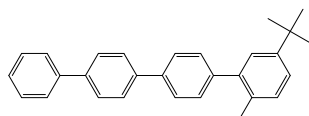


Formula	: C <sub>28</sub> H <sub>28</sub>	Dye Laser Characteristics				
MW	: 376.51 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 114932-35-3	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 290 nm(p-dioxane)	Pump source (nm)				
Fluorescence (λ-max)	: 364 nm(p-dioxane)	XeCl(308)	359	346-377	0.23	p-dioxane

Reference : 1.Appl. Phys. B32, 9(1983)

LT-D1205C

2,2"-dimethyl-1,1':4'1"-terphenyl

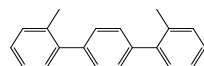


Formula	: C <sub>20</sub> H <sub>18</sub>	Dye Laser Characteristics				
MW	: 258.36 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 53092-64-1	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 266 nm(cyclohexane)	Pump source (nm)				
Fluorescence (λ-max)	: 336 nm(cyclohexane)	KrF(248)	332	311-355	0.25	Cyclohexane

Reference : 1.Appl. Phys. 20, 283(1979)

LT-D1206C


2",3,3'""-tetramethyl-1,1':4'1":4"1""-quaterphenyl

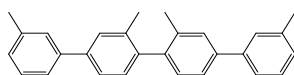

Formula	: C <sub>28</sub> H <sub>28</sub>	Dye Laser Characteristics				
MW	: 362.52 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 4575-13-7	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 266 nm(cyclohexane)	Pump source (nm)				
Fluorescence (λ-max)	: 350 nm(cyclohexane)	XeCl(308)	354	343-365	0.73	p-dioxane
		Nd:YAG(266)	350	338-361	0.73	cyclohexane

LT-D1209C


4,4"-(1,2-ethenediyl)bis-1,1'-biphenyl

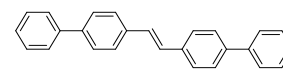

Formula	: C <sub>26</sub> H <sub>20</sub>	Dye Laser Characteristics				
MW	: 332.45 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 2039-68-1	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 341 nm(benzene)	Pump source (nm)				
Fluorescence (λ-max)	: 405 nm(benzene)	XeCl(308)	406	397-414	0.33	p-dioxane
		N2(337)	406	383-418	1	p-dioxane

LT-D1210C


2,2"-(1,1'-biphenyl)-4,4'-diyl-di-2,1-ethenediyl)bis-benzenesulfonic acid disodium salt

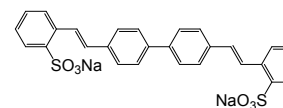

Formula	: C <sub>28</sub> H <sub>20</sub> O <sub>6</sub> S <sub>2</sub> Na <sub>2</sub>	Dye Laser Characteristics				
MW	: 562.56 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 27344-41-8	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 349 nm(methanol)	Pump source (nm)				
Fluorescence (λ-max)	: 425 nm(ethanol)	XeCl(308)	425	412-435	0.56	methanol
		N2(337)	425	414-437	0.96	methanol

LT-D1211C

1,2,3,8-tetrahydro-1,2,3,3,5-pentamethyl-7H-pyrrolo[3,2-g]quinolin-7-one

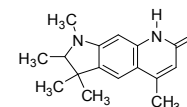


Formula	: C <sub>16</sub> H <sub>20</sub> N <sub>2</sub> O	Dye Laser Characteristics				
MW	: 256.34 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 58336-37-1	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 367 nm(ethanol)	Pump source (nm)				
Fluorescence (λ-max)	: 413 nm(ethanol)	XeCl(308)	415	399-433	0.05	methanol
		N2(337)	423	408-448	0.77	ethanol

Reference : 1.Appl. Optics, 20, 3553(1981)

LT-D1212C


4-methyl-7-(4-morpholinyl)-2H-pyran[2,3-*b*]pyridin-2-one

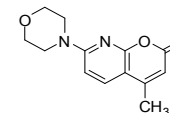

Formula	: C <sub>18</sub> H <sub>18</sub> N <sub>2</sub> O <sub>3</sub>	Dye Laser Characteristics				
MW	: 246.26 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 57980-07-1	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 357 nm(ethanol)	Pump source (nm)				
Fluorescence (λ-max)	: 428 nm(ethanol)	N2(337)	427	417-446	1.85	ethanol

LT-D1213C

7-amino-4-methyl-2H-1-benzopyran-2-one

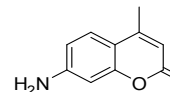


Formula	: C <sub>15</sub> H <sub>13</sub> N <sub>2</sub> O <sub>2</sub>	Dye Laser Characteristics				
MW	: 175.15 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 26093-31-2	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 354 nm(ethanol)	Pump source (nm)				
Fluorescence (λ-max)	: 428 nm(ethanol)	N2(337)	437	417-473	0.4	ethanol
		XeCl(308)	438	422-466	1.23	methanol

Reference : 1.Appl. Optics, 20, 3553(1981)

LT-D1214C

7-(ethylamino)-4-methyl-2H-1-benzopyran-2-one

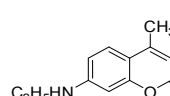


Formula	: C <sub>17</sub> H <sub>15</sub> N <sub>2</sub> O <sub>2</sub>	Dye Laser Characteristics				
MW	: 203.24 g/mole	Lasing wavelength		Concentration	Solvent	
CAS No.	: 28821-18-3	Peak(nm)	Range(nm)	(g/L)		
Absorption (λ-max)	: 363 nm(methanol)	Pump source (nm)				
Fluorescence (λ-max)	: 430 nm(ethanol)	FL	445	423-445		MeOH/H <sub>2</sub> O =1/1

Materials are used by qualified for testing and research only, there are not guaranteed in patent contention by customer use.

Materials are used by qualified for testing and research only, there are not guaranteed in patent contention by customer use.

## LT-D1215C

## 7-(ethylamino)-4,6-dimethyl-2H-1-benzopyran-2-one

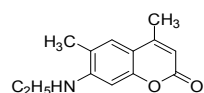


Formula : C<sub>13</sub>H<sub>15</sub>NO<sub>2</sub>  
MW : 217 g/mole  
CAS No. : 26078-25-1  
Absorption (λ-max) : 366 nm(ethanol)  
Fluorescence (λ-max) : 435 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	450	427-488	0.04	methanol
XeCl(308)	451	430-481	1.6	methanol
Nd:YAG(355)	452	436-468	0.065	methanol

Reference : 1.Appl. Optics, 13(10), 2317(1974)

## LT-D1216C

## 7-hydroxy-4-methylcoumarin

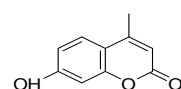


Formula : C<sub>10</sub>H<sub>8</sub>O<sub>3</sub>  
MW : 176.17 g/mole  
CAS No. : 90-33-5  
Absorption (λ-max) : 372 nm(ethanol)  
Fluorescence (λ-max) : 445 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	454		0.15	ethanol

## LT-D1217C

## 7-(diethylamino)-4-methyl-2H-1-benzopyran-2-one

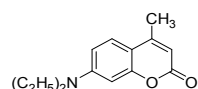


Formula : C<sub>14</sub>H<sub>17</sub>NO<sub>2</sub>  
MW : 231.3 g/mole  
CAS No. : 91-44-1  
Absorption (λ-max) : 373 nm(ethanol)  
Fluorescence (λ-max) : 446 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	457	450-484	0.035	ethanol
XeCl(308)	457	440-478	0.23	ethanol
Nd:YAG(355)	460	442-480	0.25	ethanol

Reference : 1.Appl. Optics, 24(7), 937(1985)

## LT-D1218C

## 7-(dimethylamino)-4-methyl-2H-1-benzopyran-2-one

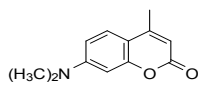


Formula : C<sub>13</sub>H<sub>13</sub>NO<sub>2</sub>  
MW : 203.24 g/mole  
CAS No. : 87-01-4  
Absorption (λ-max) : 366 nm(ethanol)  
Fluorescence (λ-max) : 453 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL		448-466	0.045	ethanol

## LT-D1219C

## 7-(diethylamino)-2H-1-benzopyran-2-one

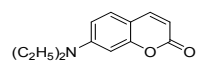


Formula : C<sub>13</sub>H<sub>15</sub>NO<sub>2</sub>  
MW : 217.26 g/mole  
CAS No. : 20571-42-0  
Absorption (λ-max) : 368 nm  
Fluorescence (λ-max) : 465 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	474	462-487	0.017	methanol
XeCl(308)	470	449-505	0.54	methanol
Nd:YAG(355)	462	450-485	0.22	ethanol

## LT-D1220C

## 2,3,6,7,10,11-hexahydro-1H,5H-cyclopenta[3,4][1]benzopyrano[6,7,8-ij]quinolizin-12(9H)-one

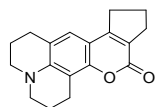


Formula : C<sub>18</sub>H<sub>18</sub>NO<sub>2</sub>  
MW : 281.36 g/mole  
CAS No. : 41175-45-5  
Absorption (λ-max) : 386 nm(ethanol)  
Fluorescence (λ-max) : 465 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	478			ethanol

## LT-D1221C

## 2,3,6,7-tetrahydro-9-methyl-1H,5H,11H-[1]benzopyrano-[6,7,8-ij]quinolizin-11-one

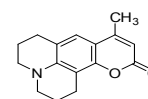


Formula : C<sub>18</sub>H<sub>17</sub>NO<sub>2</sub>  
MW : 255.32 g/mole  
CAS No. : 41267-76-9  
Absorption (λ-max) : 390 nm(ethanol)  
Fluorescence (λ-max) : 468 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	474	456-505	0.04	ethanol
XeCl(308)	476	461-506	0.77	methanol

Reference : 1.NEW Durham, NH03855

## LT-D1222C

## 7-(diethylamino)-4-(trifluoromethyl)-2H-1-benzopyran-2-one

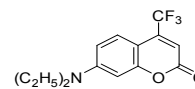


Formula : C<sub>14</sub>H<sub>14</sub>F<sub>3</sub>NO<sub>2</sub>  
MW : 285.26 g/mole  
CAS No. : 41934-47-8  
Absorption(λ-max) : 404 nm(methanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	481	475-489	0.04	p-Dioxane
XeCl(308)	485	461-515	3.3	p-Dioxane

## LT-D1223C

## 7-(dimethylamino)-4-(trifluoromethyl)-2H-1-benzopyran-2-one

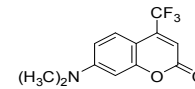


Formula : C<sub>12</sub>H<sub>10</sub>F<sub>3</sub>NO<sub>2</sub>  
MW : 257.21 g/mole  
CAS No. : 53518-14-2  
Absorption (λ-max) : 397 nm(ethanol)  
Fluorescence (λ-max) : 511 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	479			p-Dioxane
XeF(351)	519	494-566	2.6	ethanol

## LT-D1224C

## 6,7,8,9-tetrahydro-6,8,9-trimethyl-4-(trifluoromethyl)-2H-pyrano[2,3-b][1,8]naphthyridin-2-one

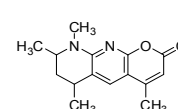


Formula : C<sub>18</sub>H<sub>15</sub>F<sub>3</sub>N<sub>2</sub>O<sub>2</sub>  
MW : 312.29 g/mole  
CAS No. : 62377-37-1  
Absorption (λ-max) : 392 nm(ethanol)  
Fluorescence (λ-max) : 480 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
XeCl(308)	489	471-535	1.66	methanol

## LT-D1225C

## 7-amino-4-(trifluoromethyl)-2H-1-benzopyran-2-one

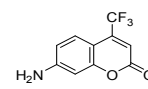


Formula : C<sub>10</sub>H<sub>7</sub>F<sub>3</sub>NO<sub>2</sub>  
MW : 229.16 g/mole  
CAS No. : 53518-15-3  
Absorption (λ-max) : 381 nm(ethanol)  
Fluorescence (λ-max) : 482 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	490	467-510	0.05	methanol

Reference : 1.Appl. Optics, 17(20), 3268(1978)

## LT-D1226C

## 2,3,6,7-tetrahydro-1H,5H,11H-[1]benzopyrano[6,7,8-ij]quinolizin-11-one

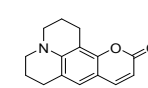

Formula : C<sub>18</sub>H<sub>15</sub>NO<sub>2</sub>  
MW : 241.37 g/mole  
CAS No. : 58336-35-9  
Absorption (λ-max) : 396 nm(methanol)  
Fluorescence (λ-max) : 475 nm(methanol)


Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	480	460-523	0.024	ethanol

Materials are used by qualified for testing and research only, there are not guaranteed in patent contention by customer use.

Materials are used by qualified for testing and research only, there are not guaranteed in patent contention by customer use.

## LT-D1227C

## 7-(ethylamino)-4-(trifluoromethyl)-2H-1-benzopyran-2-one

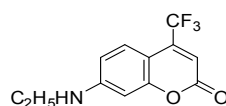


Formula : C<sub>12</sub>H<sub>10</sub>F<sub>3</sub>NO<sub>2</sub>  
MW : 257.21 g/mole  
CAS No. : 52840-38-7  
Absorption (λ-max) : 392 nm(methanol)  
Fluorescence (λ-max) : 495 nm(methanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
XeCl(308)	506	478-568	2.1	methanol
Nd:YAG(355)	500	494-504	0.26	ethanol

## LT-D1228C

## 7-(ethylamino)-6-methyl-4-(trifluoromethyl)-2H-1-benzopyran-2-one

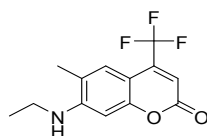


Formula : C<sub>13</sub>H<sub>12</sub>F<sub>3</sub>NO<sub>2</sub>  
MW : 271.24 g/mole  
CAS No. : 55804-70-1  
Absorption (λ-max) : 395 nm(methanol)  
Fluorescence (λ-max) : 490 nm(methanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
XeCl(308)	500	480-537	1.9	methanol
Nd:YAG(355)	500	480-541	0.4	ethanol

## LT-D1229C

## 2,3,6,7-tetrahydro-11-oxo-1H,5H,11H-[1]benzopyrano[6,7,8-ij]quinolizine-10-carboxylic acid, ethyl ester

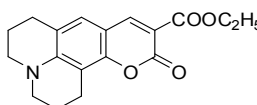


Formula : C<sub>18</sub>H<sub>19</sub>NO<sub>2</sub>  
MW : 313.34 g/mole  
CAS No. : 55804-66-5  
Absorption (λ-max) : 437 nm(methanol)  
Fluorescence (λ-max) : 479 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	505	496-519		MeOH/H <sub>2</sub> O =1/1

## LT-D1230C

## 2,3,6,7-tetrahydro-10-(3-pyridinyl)-1H,5H,11H-[1]benzopyrano[6,7,8-ij]quinolizine-11-one

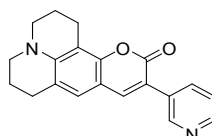


Formula : C<sub>20</sub>H<sub>18</sub>N<sub>2</sub>O<sub>2</sub>  
MW : 318 g/mole  
CAS No. : 87349-92-6  
Absorption (λ-max) : 426 nm(ethanol)  
Fluorescence (λ-max) : 496 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	511	505-518		ethanol

Reference : 1.Optic Commun., 47(1) 57(1983)

## LT-D1231C


## 3-(2-benzothiazolyl)-7-(diethylamino)-2H-1-benzopyran-2-one

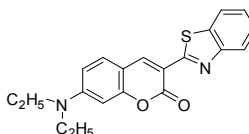

Formula : C<sub>20</sub>H<sub>18</sub>N<sub>2</sub>O<sub>2</sub>S  
MW : 350.44 g/mole  
CAS No. : 38215-36-0  
Absorption (λ-max) : 460 nm(ethanol)  
Fluorescence (λ-max) : 507 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	531	511-556	0.035	methanol methanol

## LT-D1232C

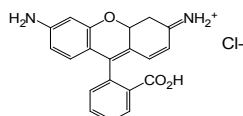
2-(6-amino-3-imino-3H-xanthen-9-yl)-benzoic acid, monohydrochloride;  
2-(6-amino-3-imino-3H-xanthen-9-yl)-benzoic acid

Formula : C<sub>20</sub>H<sub>14</sub>N<sub>2</sub>O<sub>3</sub> · HCl  
MW : 366.8 g/mole  
CAS No. : 13558-31-1  
Absorption (λ-max) : 508 nm(methanol)  
Fluorescence (λ-max) : 532 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	554	530-580	0.18	methanol
XeCl(308)	555	544-580	1.1	methanol

## LT-D1233C

## 2-[4-[4-(dimethylamino)phenyl]-1,3-butadienyl]-1,3,3-trimethyl-3H-indolium perchlorate

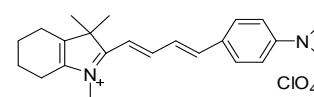


Formula : C<sub>23</sub>H<sub>27</sub>N<sub>2</sub> · ClO<sub>4</sub><sup>-</sup>  
MW : 430.93 g/mole  
CAS No. : 76433-27-7  
Absorption (λ-max) : 605 nm(methanol)  
Fluorescence (λ-max) : 670 nm(methanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
Nd:YAG(532)	716	692-743		methanol

## LT-D1234C

## 2-[4-[4-(dimethylamino)phenyl]-1,3-butadienyl]-3-ethyl-naphtho[2,1-c]thiazolium perchlorate

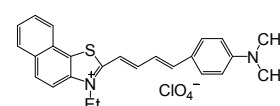


Formula : C<sub>25</sub>H<sub>25</sub>N<sub>2</sub>S · ClO<sub>4</sub><sup>-</sup>  
MW : 484 g/mole  
CAS No. : 89872-07-1  
Absorption (λ-max) : 573 nm(methanol)  
Fluorescence (λ-max) : 706 nm(methanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
Nd:YAG(532)	718	700-743	0.15	methanol

## LT-D1235C

## 4-[4-[4-(dimethylamino)phenyl]-1,3-butadienyl]-1-ethyl quinolinium perchlorate

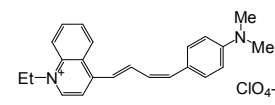


Formula : C<sub>23</sub>H<sub>25</sub>N<sub>2</sub> · ClO<sub>4</sub><sup>-</sup>  
MW : 428.9 g/mole  
CAS No. : 92479-59-9  
Absorption (λ-max) : 558 nm(methanol)  
Fluorescence (λ-max) : 766 nm(methanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
Nd:YAG(532)	785	758-826	0.14	ethanol

## LT-D1236C

## 2-(6-(9-(2,3,6,7-tetrahydro-1H,5H-benzo(i,j)quinolizinium))-2,4-neopentylene-1,3,5-hexatrienyl)-3-methylbenzothiazolium perchlorate

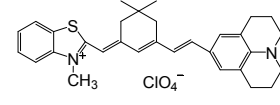


Formula : C<sub>31</sub>H<sub>35</sub>N<sub>2</sub>S · ClO<sub>4</sub><sup>-</sup>  
MW : 567.15 g/mole  
CAS No. : 385437-35-4

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
Nd:YAG(532)	880	856-918	0.57	propylene carbonate
XeCl(308)	881	856-919	1.14	DMSO

## LT-D1237C

## 2-[5-(1,3-dihydro-1,3,3-trimethyl-2H-indol-2-ylidene)-1,3-pentadienyl]-1,3,3-trimethyl-3H-indolium iodide or (1,1',3,3,3',3'-hexamethyl-2,2'-indodicarbocyanine iodide)

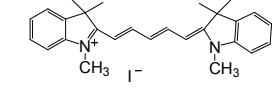


Formula : C<sub>27</sub>H<sub>31</sub>N<sub>2</sub> · I<sup>-</sup>  
MW : 510.46 g/mole  
CAS No. : 36536-22-8  
Absorption (λ-max) : 643 nm(ethanol)  
Fluorescence (λ-max) : 670 nm(ethanol)

Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
Nd:YAG(585)	675		0.1	methanol

## LT-D1238C

## 3-ethyl-2-[7-(3-ethyl-2(3H)-benzoxazolylidene)-1,3,5-heptatrienyl]-benzoxazolium iodide

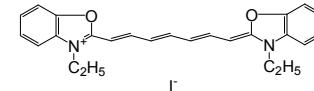

Formula : C<sub>25</sub>H<sub>25</sub>N<sub>2</sub>O<sub>2</sub> · I<sup>-</sup>  
MW : 512.39 g/mole  
CAS No. : 15185-43-0  
Absorption (λ-max) : 686 nm(ethanol)  
Fluorescence (λ-max) : 726 nm(ethanol)


Pump source (nm)	Lasing wavelength		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
XeCl(308)	798	777-800	0.77	DMSO
N2(337)	777	762-802	0.61	DMSO

Materials are used by qualified for testing and research only, there are not guaranteed in patent contention by customer use.

Materials are used by qualified for testing and research only, there are not guaranteed in patent contention by customer use.

LT-D1240C

2-[7-(1,3-dihydro-1,3,3-trimethyl-2H-indol-2-ylidene)-1,3,5-heptatrienyl]-1,3,3-trimethyl-3H-indolium iodide

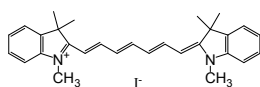


Formula :  $C_{29}H_{33}N_2 \cdot I^-$   
MW : 536.5 g/mole  
CAS No. : 19764-96-6  
Absorption ( $\lambda$ -max) : 751 nm(DMSO)  
Fluorescence ( $\lambda$ -max) : 792 nm(DMSO)

Pump source (nm)	Dye Laser Characteristics		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
XeCl(308)	826	812-838	0.75	methanol
XeCl(308)	851	822-887	1.23	DMSO
N2(337)	846	828-891	1	DMSO

LT-D1241C

2-[7-(1,3-dihydro-1,3,3-trimethyl-2H-indol-2-ylidene)-1,3,5-heptatrienyl]-1,3,3-trimethyl-3H-indolium perchlorate

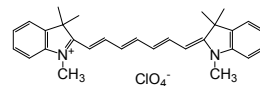


Formula :  $C_{29}H_{33}N_2 \cdot ClO_4^-$   
MW : 509.04 g/mole  
CAS No. : 16595-48-5  
Absorption ( $\lambda$ -max) : 750 nm(DMSO)  
Fluorescence ( $\lambda$ -max) : 790 nm(DMSO)

Pump source (nm)	Dye Laser Characteristics		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
XeCl(308)	826	812-838	0.75	methanol
XeCl(308)	851	822-887	1.23	DMSO
N2(337)	846	828-890	1	DMSO

LT-D1242C

3-ethyl-2-[7-(3-ethyl-2(3H)-benzothiazolylidene)-1,3,5-heptatrienyl]-benzothiazolium iodide

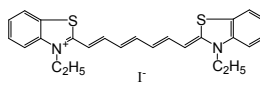


Formula :  $C_{25}H_{25}N_2S_2 \cdot I^-$   
MW : 544.51 g/mole  
CAS No. : 3071-70-3  
Absorption ( $\lambda$ -max) : 761 nm(ethanol)  
Fluorescence ( $\lambda$ -max) : 813 nm(ethanol)

Pump source (nm)	Dye Laser Characteristics		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
XeCl(308)	849	828-883	0.42	DMSO
XeCl(308)	879	836-925	0.38	DMSO
N2(337)	855	840-877	0.65	DMSO

LT-D1243C

2-[7-(1,3-dihydro-1,1-dimethyl-3-(4-sulfobutyl)-2H-benz[e]indol-2-ylidene)-1,3,5-heptatrienyl]-1,1-dimethyl-3-(4-sulfobutyl)-1H-benz[e]indolium hydroxide, inner salt, sodium salt

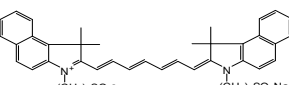


Formula :  $C_{33}H_{48}N_2O_6S_2 \cdot Na$   
MW : 774.97 g/mole  
CAS No. : 3599-32-4  
Absorption ( $\lambda$ -max) : 795 nm(DMSO)  
Fluorescence ( $\lambda$ -max) : 832 nm(DMSO)

Pump source (nm)	Dye Laser Characteristics		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
Nd:YAG(532)	863	845-905	0.08	DMSO
N2(337)	887	872-935	1	DMSO

LT-D1244C

2-[2-(2-(diphenylamino)-3-[(1-ethylnaphtho[1,2-d]thiazol-2(H)-ylidene)ethylidene]-1-cyclopenten-1-yl)ethenyl]-1-ethyl-naphtho[1,2-d]thiazolium perchlorate

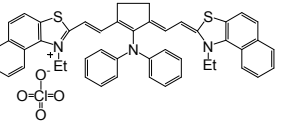


Formula :  $C_{47}H_{40}N_2S_2 \cdot ClO_4^-$   
MW : 810.42 g/mole  
CAS No. : 54849-65-9  
Fluorescence ( $\lambda$ -max) : 839 nm(methanol)

Pump source (nm)	Dye Laser Characteristics		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	972			

LT-D1245C

5-chloro-2-[2-[3-[(5-chloro-3-ethyl-2(3H)-benzothiazol-ylidene)ethylidene]-2-(diphenylamino)-1-cyclopenten-1-yl]ethenyl]-3-ethyl benzothiazolium perchlorate

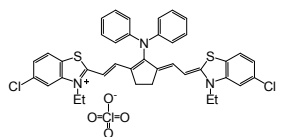


Formula :  $C_{39}H_{34}Cl_2N_2S_2 \cdot ClO_4^-$   
MW : 779.2 g/mole  
CAS No. : 53655-17-7  
Absorption ( $\lambda$ -max) : 822 nm(DMSO)  
Fluorescence ( $\lambda$ -max) : 881 nm(DMSO)

Pump source (nm)	Dye Laser Characteristics		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	950			
XeCl(308)	880	868-900	0.25	DMF
N2(337)	912	896-939	1.17	DMSO

LT-D1246C

3-ethyl-2-[5-(3-ethyl-2(3H)-benzoxazolylidene)-1,3-pentadienyl]-benzoxazolium iodide

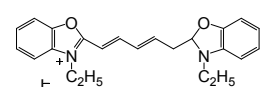


Formula :  $C_{23}H_{23}N_2O_2 \cdot I^-$   
MW : 486.35 g/mole  
CAS No. : 14806-50-9  
Absorption ( $\lambda$ -max) : 582 nm(ethanol)  
Fluorescence ( $\lambda$ -max) : 610 nm(ethanol)

Pump source (nm)	Dye Laser Characteristics		Concentration (g/L)	Solvent
	Peak(nm)	Range(nm)		
FL	662			DMSO

LT-D1248C


2-[2-[2-chloro-3-[(1,3-dihydro-1,3,3-trimethyl-2H-indol-2-ylidene)-ethylidene]-1-cyclohexen-1-yl]ethenyl]-1,3,3-trimethylindolium iodide

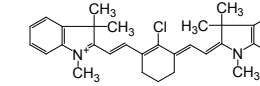

Formula :  $C_{32}H_{36}ClN_2 \cdot I^-$   
MW : 611.02 g/mole  
Absorption ( $\lambda$ -max) : 775 nm(methanol)  
Fluorescence ( $\lambda$ -max) : 817 nm(ethanol)

LT-D1249C


2-[2-[2-chloro-3-[(1,3-dihydro-1,3,3-trimethyl-2H-indol-2-ylidene)-ethylidene]-1-cyclohexen-1-yl]ethenyl]-1,3,3-trimethyl-1H-indolium perchlorate

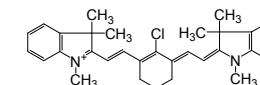

Formula :  $C_{32}H_{36}ClN_2 \cdot ClO_4^-$   
MW : 583.56 g/mole  
Absorption ( $\lambda$ -max) : 776 nm(methanol)

LT-D1250C


2-[2-[2-chloro-3-[(1,3-dihydro-3,3-dimethyl-1-propyl-2H-indol-2-ylidene)-ethylidene]-1-cyclohexen-1-yl]ethenyl]-3,3-dimethyl-1-propylindolium iodide

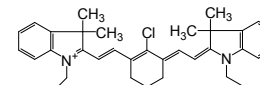

Formula :  $C_{38}H_{44}ClN_2 \cdot I^-$   
MW : 667.12 g/mole  
Absorption ( $\lambda$ -max) : 776 nm(methanol)  
Fluorescence ( $\lambda$ -max) : 817 nm

LT-D1251C


2-[2-[2-chloro-3-[(1,3-dihydro-3,3-dimethyl-1-propyl-2H-indol-2-ylidene)-ethylidene]-1-cyclohexen-1-yl]ethenyl]-3,3-dimethyl-1-propyl-1H-indolium perchlorate

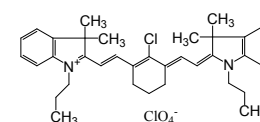

Formula :  $C_{38}H_{44}ClN_2 \cdot ClO_4^-$   
MW : 639.66 g/mole  
Absorption ( $\lambda$ -max) : 778 nm(methanol)

LT-D1253C

2-[2-[2-chloro-3-[(1,3-dihydro-3,3-dimethyl-1-ethyl-2H-benz[e]indol-2-ylidene)-ethylidene]-1-cyclohexen-1-yl]ethenyl]-3,3-dimethyl-1-ethyl-1H-benz[e]indolium iodide

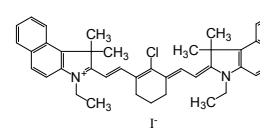

Formula :  $C_{42}H_{44}ClN_2 \cdot I^-$   
MW : 739.18 g/mole  
Absorption ( $\lambda$ -max) : 815 nm(methanol)

Materials are used by qualified for testing and research only, there are not guaranteed in patent contention by customer use.

Materials are used by qualified for testing and research only, there are not guaranteed in patent contention by customer use.